

Trade fairs cancelled. And now? LCTech presents its product news.

BE PART OF IT



Login 

German  

Print 

Search

LCTech-en

Pickering

Maintenance Onyx PCX

Maintenance Onyx PCX



Servicing of Your Onyx PCX

Keep your Onyx PCX in good condition and minimize downtime by performing annual maintenance.

Below you can find a general maintenance schedule. Your individual maintenance schedule will depend on usage. If you use the Onyx PCX very intensely, you may want to perform the preventative maintenance more often.

Refer to the maintenance section of the Pinnacle manual for details on performing the maintenance. The manual is ready for download on the

sidebar.

For professional maintenance complete maintenance and tubing kits are available.

Routine maintenance

- Check for leaks every day
- Replace flush and piston wash reservoirs every week
- Record operating pressures in a daily log book

Precaution is cheaper than repair

- Replace valve and pump seals every 12 months
- Replace the 10 µm reagent frits every 3 months
- Change the tubing's every 24 months

Please contact us

+49 8082  

2717-0info@LCTech.de

Recommendations for Maintenance

Onyx PCX

[Recommendations_for_Maintenanc
e_Onyx_PCX.pdf \(pdf | 297 KB \)](#)

Spare parts for Onyx PCX

[Overview: Spare parts for Onyx
PCX \(external link\)](#)

Manual

[Manual Onyx PCX \(pdf | 3 MB \)](#)

© 2020 LCTech GmbH

[Disclaimer](#) |
[Data Security Statement](#) |

[Terms of Service](#)

[Sitemap](#)